




D-NET WITH RETRACTABLE POLE

Code 0168

1. Slide the shaft of the D-net into the inner pole of the retractable pole.
2. Align the holes on the D-net shaft with the holes on the retractable pole.
3. Place one curved washer on the screw.
4. Thread the screw through the holes in the D-net shaft and the pole.
5. Place the second curved washer over the other end of the screw as shown.
6. For the best results, use a power screwdriver to drive the screw through the second hole and tighten the washer to the pole.
7. To extend the pole, unscrew the large white collar. Extend the pole to the desired length. Slide the large white collar along the pole to meet the small white collar. Screw the large white collar onto the small white collar to lock the pole at the desired length. **Do not over extend pole, as inner pole will come out of outer pole.**

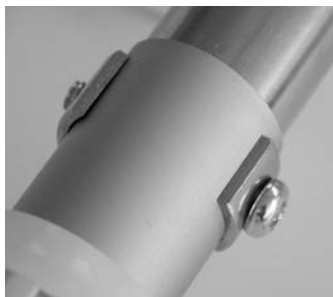

LaMOTTE COMPANY

PO Box 329 • Chestertown • Maryland • 21620 • USA
800-344-3100 • 410-778-3100 (Outside U.S.A.) • Fax 410-778-6394
Visit us on the web at www.lamotte.com